

Evaluating product spreading characteristics on a rotational rheometer using the power law model



RHEOLOGY AND VISCOSITY

Introduction

The rheological properties of a product can influence how it is visually and texturally perceived by a consumer and how it is likely to behave during product use. For example very shear thinning materials will be highly responsive to changes in applied stress while Newtonian materials will show much lesser dependence. Such a response is important when considering the ease of spreading or 'spread-ability'.

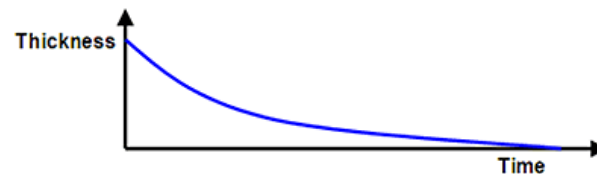


Figure 1: Diagram indicating how product layer thickness changes during spreading

The process of spreading causes a consequent reduction in the layer thickness as it is distributed over a wider surface area as shown in Figure 1. Since shear rate is equal to the applied velocity divided by the layer thickness, spreading can therefore not be attributed to a single shear rate.

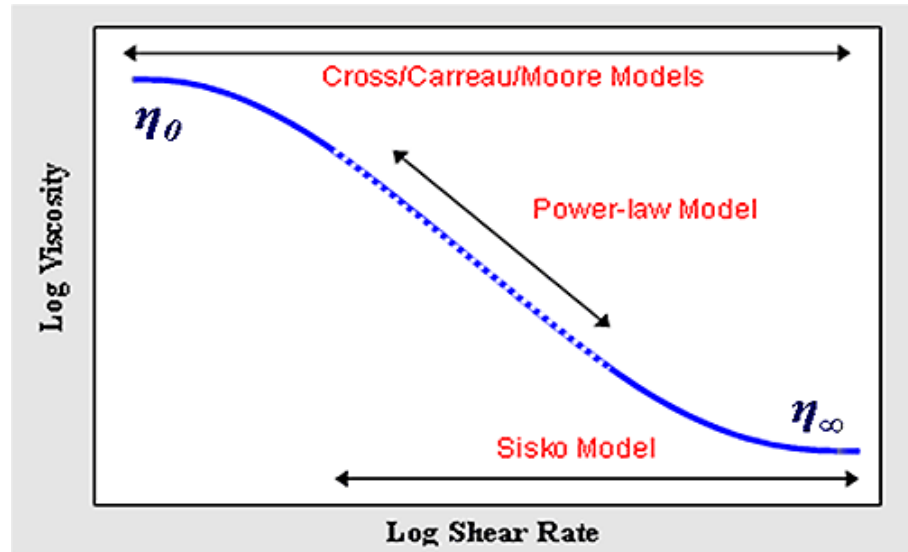


Figure 2: Diagram showing an ideal flow curve and the relevant models for describing its shape

A better way of assessing spread-ability is by characterizing the change in viscosity over a range of shear rates as shown in Figure 2. The region of interest is the shear thinning or power law region since this describes how easily the material structure breaks down with applied shear. This region appears linear on a log-log plot of viscosity vs. shear rate with a constant gradient, but shows power law dependence when plotted on a linear scale.

Mathematically this region of the flow curve can be described using the Power Law or Ostwald de Waele Model given by Equation 1.

$$\sigma = k\dot{\gamma}^n$$

k is the consistency

n is the power law index

σ is the shear rate

$\dot{\gamma}$ is the shear rate

Consistency has the units of Pas^n but is numerically equal to the viscosity measured at 1s^{-1} . The power law index ranges from 0 for very shear thinning materials to 1 for Newtonian materials.

The lower the stress input required the easier the material should be to spread. A lower value of k means lower viscosity and hence lower stress input, while a lower value of n indicates greater shear thinning which translates to a smaller stress increase with increasing shear rate.

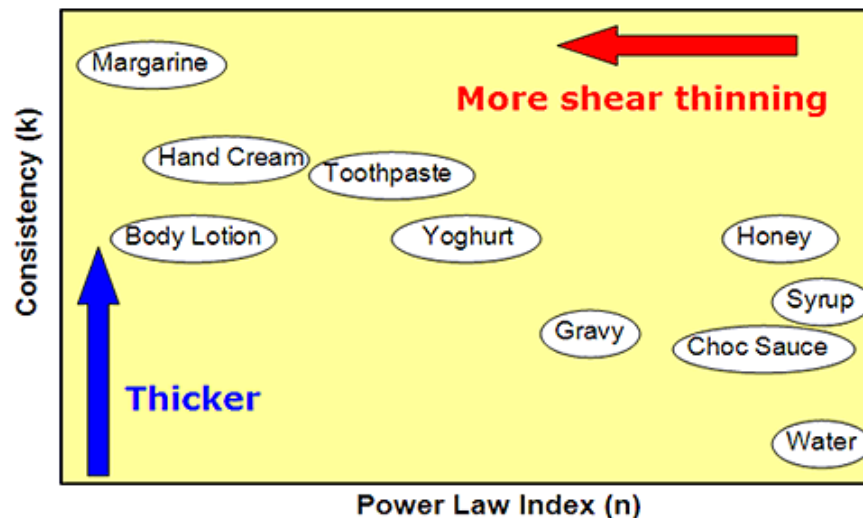


Figure 3: Diagram showing how different products may fit on a plot of k versus n

This information can be represented on a chart similar to that shown in Figure 3. Materials with low k values and/or low n values should be easiest to spread.

Experimental

- The spread-ability of a number of consumer products was evaluated by performing a shear rate ramp test and analyzing the resultant curve using a power law model.
- Rotational rheometer measurements were made using a Kinexus rotational rheometer with a Peltier plate cartridge and a roughened parallel plate measuring system¹, and using standard pre-configured sequences in the rSpace software.
- A standard loading sequence was used to ensure that both samples were subject to a consistent and controllable loading protocol.
- All rheology measurements were performed at 25°C.
- A flow curve was generated using a shear rate ramp test and a power law model fitted to this curve.

Results and Discussion

Figure 4 shows the viscosity-shear rate curve for a number of commercial products and their corresponding fitting parameters, with graphical presentation of the latter in Figure 5.

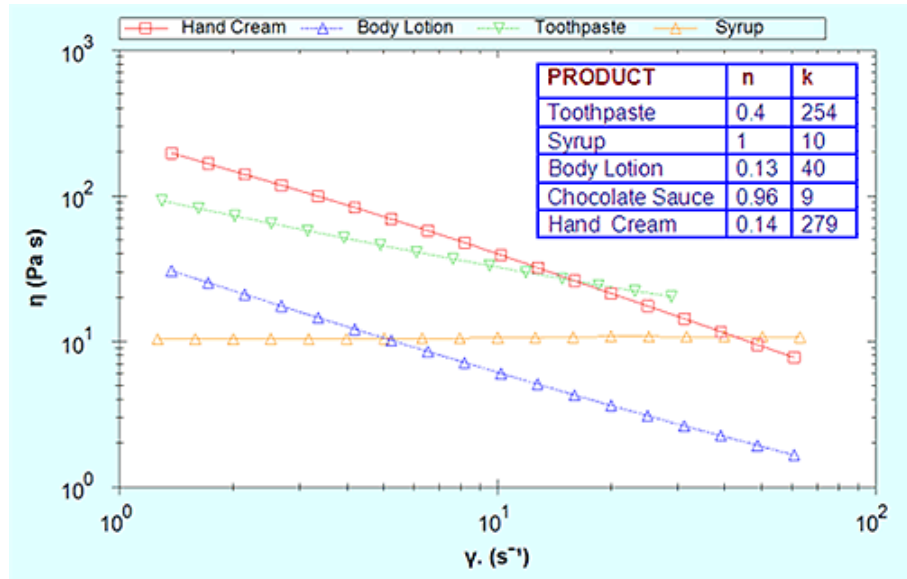


Figure 4: Flow curves and model fitting parameters for various samples

Although toothpaste and hand cream have similar k values, hand cream has a much lower n value making it more shear thinning and easier to spread. Conversely syrup and chocolate sauce have much lower k values but are not shear thinning hence they appear thick and sticky during application. The body lotion has both a relatively low k and n value making it much easier to apply.

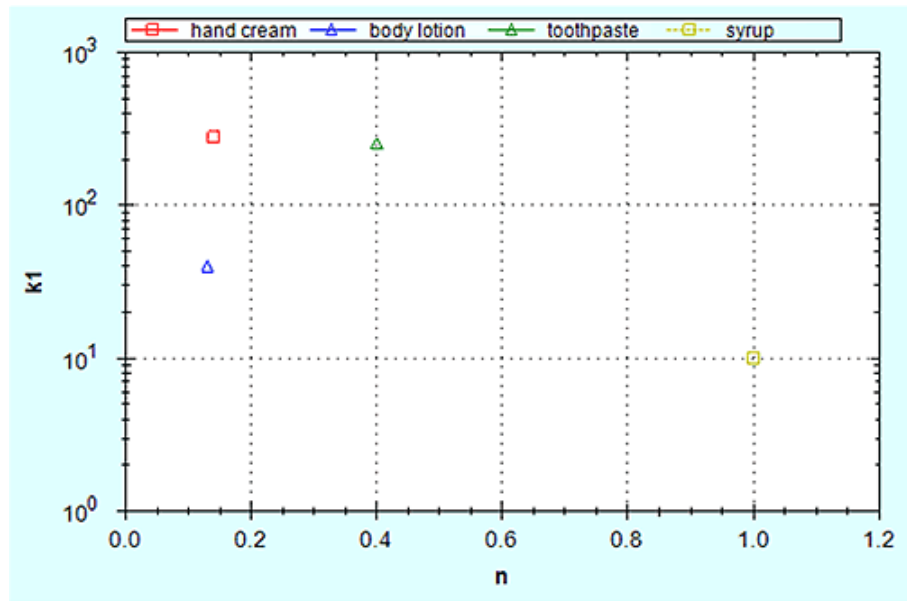


Figure 5: Model fitting parameters k and n plotted against one another

To quantitatively compare the stress requirement for the spreading of a hand cream and syrup respectively at equivalent shear rates the values of n and k can be substituted into Equation 1. Considering a single shear rate of 1s^{-1} which may equivocate with a thicker product layer then the stress required to maintain flow at this shear rate is 279Pa for the hand cream and 10Pa for syrup ($\sigma = k$ at 1s^{-1}). At a shear rate of 1000s^{-1} which would relate to a thinner layer of material resulting from the

spreading process the stress requirement increases to 734Pa for the hand cream and 10,000Pa for the syrup. This highlights importance of non-Newtonian behaviour in the spreading process.

Conclusion

A shear rate ramp test with a power law model fit was used to characterize the spreadability of various commercial products using the power law fitting parameters k and n . Low values of k and n indicate lower viscosity and a greater degree of shear thinning respectively which will contribute to easier spreading.

Notes

1. Note that testing is recommended to be undertaken with cone and plate or parallel plate geometry – with the latter being preferred for dispersions and emulsions with large particle sizes. Such material types may also require the use of serrated or roughened geometries to avoid artefacts relating to slippage at the geometry surface.



Malvern Instruments Limited
Groewood Road, Malvern,
Worcestershire, UK. WR14
1XZ

Tel: +44 1684 892456
Fax: +44 1684 892789
www.malvern.com

Malvern Instruments is part of Spectris plc, the Precision Instrumentation and Controls Company.

Spectris and the Spectris logo are Trade Marks of Spectris plc.

spectris

All information supplied within is correct at time of publication.

Malvern Instruments pursues a policy of continual improvement due to technical development. We therefore reserve the right to deviate from information, descriptions, and specifications in this publication without notice. Malvern Instruments shall not be liable for errors contained herein or for incidental or consequential damages in connection with the furnishing, performance or use of this material.

Malvern Instruments owns the following registered trademarks: Bohlin, FIPA, Insitex, ISYS, Kinexus, Malvern, Malvern 'Hills' logo, Mastersizer, MicroCal, Morphologi, Rosand, 'SEC-MALS', Viscosizer, Viscotek, Viscogel and Zetasizer.